



# PFAS on Force C18 by Direct Injection (EPA Method 8327)

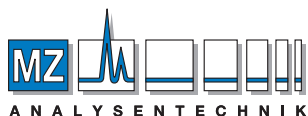

AUTHORIZED DISTRIBUTOR

MZ-Analysentechnik GmbH  
 Barcelona-Allee 17 • D-55129 Mainz  
 Tel +49 6131 880 96-0  
 Fax +49 6131 880 96-20  
 e-mail: info@mz-at.de  
 www.mz-at.de

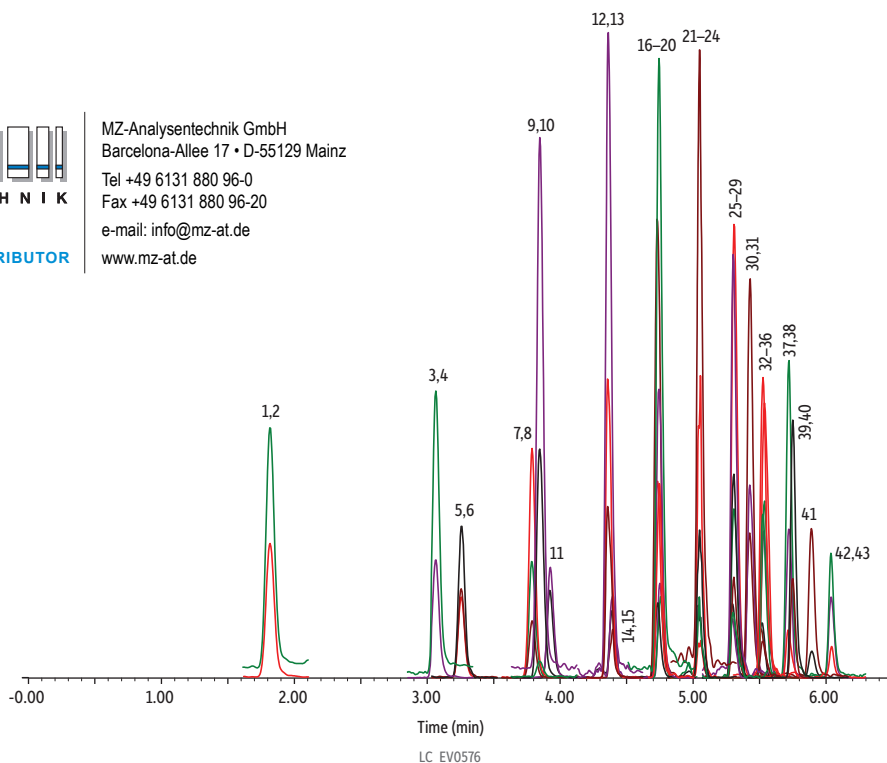

Peaks	tr (min)	Conc. (ng/L)	Precursor Ion	Product Ion	Product Ion	Column	Force C18 (cat.# 9634252)
1. Perfluorobutanoic acid (PFBA)	1.82	80	231.1	169.1	-	Dimensions:	50 mm x 2.1 mm ID
2. Perfluoro- <i>n</i> -[ <sup>13</sup> C <sub>4</sub> ]butanoic acid (M4PFBA)	1.82	40	217.0	172.0	-	Particle Size:	1.8 μm
3. Perfluoropentanoic acid (PFPeA)	3.07	80	263.0	219.0	-	Pore Size:	100 Å
4. Perfluoro- <i>n</i> -[ <sup>13</sup> C <sub>5</sub> ]pentanoic acid (M5PFPeA)	3.07	40	268.0	223.0	-	Temp.:	40 °C
5. Perfluorobutane sulfonic acid (PFBS)	3.25	80	299.1	80.0	99.0	Sample	
6. Perfluoro-1-[2,3,4- <sup>13</sup> C <sub>3</sub> ]butyl sulfonic acid (M3PFBS)	3.25	40	302.0	80.0	-	Diluent:	50:50 Water:methanol, 0.1% acetic acid
7. 1H, 1H, 2H, 2H-perfluorohexane sulfonic acid (4:2 FTS)	3.79	80	327.1	307.1	80.8	Conc.:	80 ppt (40 ppt for surrogate or isotopically labelled PFAS)
8. 1H, 1H, 2H, 2H-perfluoro-1-[1,2- <sup>13</sup> C <sub>2</sub> ]hexyl sulfonic acid (M2-4:2 FTS)	3.79	40	329.0	309.0	-	Inj. Vol.:	10 μL
9. Perfluorohexanoic acid (PFHxA)	3.84	80	313.1	269.1	119.1	Mobile Phase	
10. Perfluoro- <i>n</i> -[1,2,3,4,6- <sup>13</sup> C <sub>5</sub> ]hexanoic acid (M5PFHxA)	3.84	40	318.0	273.0	-	A:	Water, 5 mM ammonium acetate
11. Perfluoropentane sulfonic acid (PFPeS)	3.92	80	349.1	80.0	99.0	B:	Methanol
12. Perfluoroheptanoic acid (PFHpA)	4.36	80	363.2	319.1	169.1	Time (min)	Flow (mL/min)
13. Perfluoro- <i>n</i> -[1,2,3,4- <sup>13</sup> C <sub>4</sub> ]heptanoic acid (M4PFHpA)	4.36	40	367.0	322.0	-	0.00	0.4
14. Perfluorohexane sulfonic acid (PFHxS)	4.40	80	399.1	80.0	99.0	6.00	0.4
15. Perfluoro-1-[1,2,3- <sup>13</sup> C <sub>3</sub> ]hexyl sulfonic acid (M3PFHxS)	4.40	40	402.0	80.0	-	6.00	0.4
16. 1H, 1H, 2H, 2H-perfluorooctane sulfonic acid (6:2 FTS)	4.72	80	427.2	407.2	80.7	6.00	0.4
17. 1H, 1H, 2H, 2H-perfluoro-1-[1,2- <sup>13</sup> C <sub>2</sub> ]octyl sulfonic acid (M2-6:2 FTS)	4.72	40	429.0	409.0	-	8.00	0.4
18. Perfluorooctanoic acid (PFOA)	4.75	80	412.9	369.0	169.0	%A	%B
19. Perfluoro- <i>n</i> -[ <sup>13</sup> C <sub>8</sub> ]octanoic acid (M8PFOA)	4.75	40	421.0	376.0	-	0.00	80
20. Perfluoroheptane sulfonic acid (PFHpS)	4.75	80	449.2	80.0	99.0	6.00	5
21. Perfluorononanoic acid (PFNA)	5.05	80	463.2	419.2	219.1	6.00	80
22. Perfluoro- <i>n</i> -[ <sup>13</sup> C <sub>9</sub> ]nonanoic acid (M9PFNA)	5.05	40	472.0	427.0	-	8.00	80
23. Perfluorooctane sulfonic acid (PFOS)	5.05	80	499.2	80.0	99.0	Detector	MS/MS
24. Perfluoro-1-[ <sup>13</sup> C <sub>8</sub> ]octyl sulfonic acid (M8PFOS)	5.05	40	506.8	80.0	-	Ion Mode:	ESI-
25. 1H, 1H, 2H, 2H-perfluorodecane sulfonic acid (8:2 FTS)	5.29	80	527.2	507.2	80.8	Mode:	MRM
26. 1H, 1H, 2H, 2H-perfluoro-1-[1,2- <sup>13</sup> C <sub>2</sub> ]decyl sulfonic acid (M2-8:2 FTS)	5.29	40	529.0	509.0	-	Instrument	UHPLC
27. Perfluorononane sulfonic acid (PFNS)	5.30	80	549.2	80.0	99.0	Notes	A PFAS delay column (cat.# 27854) was installed before the injector.
28. Perfluorodecanoic acid (PFDA)	5.30	80	513.2	469.2	219.1		
29. Perfluoro- <i>n</i> -[1,2,3,4,5,6- <sup>13</sup> C <sub>6</sub> ]decanoic acid (M6PFDA)	5.30	40	518.9	473.9	-		
30. N-methylperfluoro-1-octanesulfonamidoacetic acid (N-MeFOSAA)	5.43	80	570.2	419.2	483.2		
31. N-methyl-d <sub>3</sub> -perfluoro-1-octanesulfonamidoacetic acid (d <sub>3</sub> -N-MeFOSAA)	5.43	40	572.9	418.9	-		
32. Perfluoroundecanoic acid (PFUnA)	5.52	80	563.2	519.2	269.1		
33. Perfluoro- <i>n</i> -[1,2,3,4,5,6,7- <sup>13</sup> C <sub>7</sub> ]undecanoic acid (M7PFUnA)	5.52	40	569.9	524.9	-		
34. Perfluorodecane sulfonic acid (PFDS)	5.53	80	599.2	80.0	99.0		
35. N-ethylperfluoro-1-octanesulfonamidoacetic acid (N-EtFOSAA)	5.54	80	584.2	419.2	483.1		
36. N-ethyl-d <sub>5</sub> -perfluoro-1-catanesulfonamidoacetic acid (d <sub>5</sub> -N-EtFOSAA)	5.54	40	589.0	419.0	-		
37. Perfluorododecanoic acid (PFDoA)	5.71	80	613.2	569.2	169.1		
38. Perfluoro- <i>n</i> -[1,2- <sup>13</sup> C <sub>2</sub> ]dodecanoic acid (M2PFDoA)	5.71	40	614.8	569.9	-		
39. Perfluorooctane sulfonamide (FOSA)	5.75	80	498.2	78.0	-		
40. Perfluoro-1-[ <sup>13</sup> C <sub>8</sub> ]octanesulfonamide (M8FOSA)	5.75	40	505.9	78.0	-		
41. Perfluorotridecanoic acid (PFTriA)	5.89	80	663.2	619.2	169.1		
42. Perfluorotetradecanoic acid (PFTreA)	6.04	80	713.2	669.2	169.1		
43. Perfluoro- <i>n</i> -[1,2- <sup>13</sup> C <sub>2</sub> ]tetradecanoic acid (M2PFTreA)	6.04	40	714.8	669.7	-		



© 2021 Restek Corporation. All rights reserved.  
 www.restek.com

Restek patents and trademarks are the property of Restek Corporation. (See www.restek.com/Patents-Trademarks for full list.) Other trademarks in Restek literature or on its website are the property of their respective owners. Restek registered trademarks are registered in the U.S. and may also be registered in other countries.