


Acrylamide (Potato Chip Extract) Stabilwax®

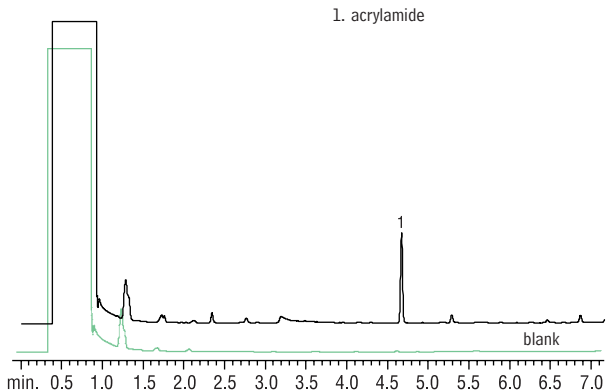

GC_FF00643

Stabilwax® 15m, 0.53mm ID, 0.50 μ m (cat.# 10637)
Sample: 1g of potato chips spiked with 100 μ g acrylamide
Inj.: 1.0 μ L, 0.5min hold
Liner: 2mm splitless with wool (cat.# 20830)
Inj. temp.: 260°C
Carrier gas: helium, constant pressure
Linear velocity: 62cm/sec. @ 100°C
Oven temp.: 100°C (hold 0.5 min.) to 200°C @ 15°C/min.
Det.: FID @ 260°C